

VICI VALCO VALVES

Clarity Control Module

ENG

Code/Rev.: M111/80C
Date: 8/24/2020

Phone: +420 251 013 400
Fax: +420 251 013 401
clarity@dataapex.com
www.dataapex.com

DataApex Ltd.
Petrzilkova 2583/13
158 00 Prague 5
The Czech Republic

Clarity[®], DataApex[®] and ▲[®] are trademarks of DataApex Ltd. Microsoft[®] and Windows[™] are trademarks of Microsoft Corporation.

DataApex reserves the right to make changes to manuals without prior notice. Updated manuals can be downloaded from www.dataapex.com.

Author: DM

Contents

1 VICI Valco Valves Control Module	1
2 Requirements	2
3 Installation Procedure	3
3.1 VICI Valco Valves Communication	3
3.2 Clarity Configuration	4
4 Using the Control Module	6
4.1 Method Setup - Valves	6
4.2 Method Setup - Event Table	8
4.3 Device Monitor	9
4.4 Valco Valve Actuator Setup	10
5 Report Setup	13
6 Troubleshooting	14

To facilitate the orientation in the **VICI Valco Valves** manual and **Clarity** chromatography station, different fonts are used throughout the manual. Meanings of these fonts are:

Instrument (blue text) marks the name of the window to which the text refers.

Open File (italics) describes the commands and names of fields in **Clarity**, parameters that can be entered into them or a window or dialog name (when you already are in the topic describing the window).

WORK1 (capitals) indicates the name of the file and/or directory.

ACTIVE (capital italics) marks the state of the station or its part.

The bold text is sometimes also used for important parts of the text and the name of the **Clarity** station. Moreover, some sections are written in format other than normal text. These sections are formatted as follows:

Note: Notifies the reader of relevant information.

Caution: Warns the user of possibly dangerous or very important information.

Marks the problem statement or trouble question.

Description: Presents more detailed information on the problem, describes its causes, etc.

Solution: Marks the response to the question, presents a procedure how to remove it.

1 VICI Valco Valves Control Module

This manual describes the setting of any **VICI Valco** valve as long as it fits on one of the following actuators: **Two-position Microelectric** actuator with RS232, **Multiposition Microelectric** actuator with RS232 or **Universal** actuator with RS232 or USB communication. The control module enables direct control of the instrument over serial line or USB port.

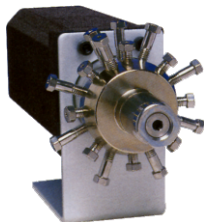


Fig 1: VICI Valco multiposition valve

Direct control means that the valve can be completely controlled from the **Clarity** environment. Instrument method, controlling the analysis conditions will be saved in the measured chromatograms.

2 Requirements

- **Clarity** Installation USB with LC Control (p/n A24) or GC Control module (p/n A23).

Note: **VICI Valco** Valves are supported since **Clarity** version **4.0.3**. In previous versions of **Clarity**, Legacy mode 2 needs to be used.

- Free serial COM (RS-232) port in the PC.

Note: Modern computers usually have only 1 (if any) serial port installed. To use more devices requiring the serial port, the **MultiCOM** adapter (p/n MC01) is available. Multiple **VICI Valco** valves still use only one serial port, but other directly controlled devices may use another serial ports.

- Serial cable (p/n SK17) provided by **DataApex** or from the valve's manufacturer (for more details see also chapter **Installation Procedure** on pg 3).

Note: Cables are not part of **Clarity** Control Module. It is strongly recommended to order required cables together with the Control Module.

- In case of USB communication, free USB port in the PC.

Note: When using USB port, USB-COM adapter drivers need to be installed. They are standard FTDI drivers available in ...\\Clarity\\BIN\\HW_DRIVERS\\FTDI folder. The configuration is same as when using RS-232 as the virtual COM port is created.

3 Installation Procedure

3.1 VICI Valco Valves Communication

The **VICI Valco** valves are controlled by serial communication. It uses special serial cable provided by the valves manufacturer (**VICI Valco** product number I-22697), but can also be ordered from **DataApex** (p/n SK17). On the computer side, it bears standard DB9F connector, on the valve's side there is special 3-pin connector. The wiring of the cable is shown on the picture:

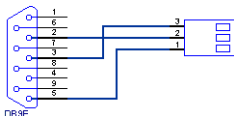


Fig 2: Serial cable for VICI Valco valves direct control

The communication parameters are:

Baud rate 9600, parity N, bits 8, stop bit 1.

3.2 Clarity Configuration

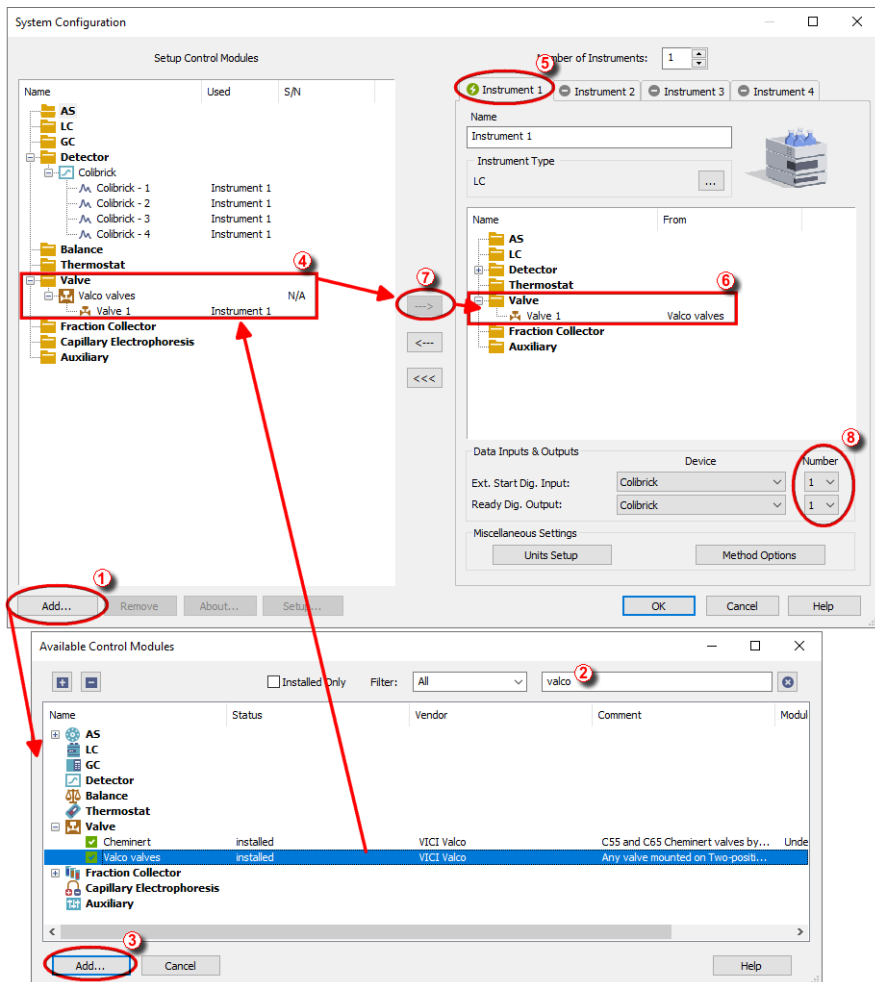



Fig 3: System Configuration

- Start the **Clarity** station by clicking on the  icon on the desktop.
- Invoke the **System Configuration** dialog accessible from the **Clarity** window using the **System - Configuration...** command.
- Press the **Add** button ① (see **Fig 3** on pg 4.) to invoke the **Available Control Modules** dialog.
- You can specify the searching filter ② to simplify the finding of the driver.

- Select the **ET, EH, EP, EHM, EMT, ECHM, ECMT** item and press the **Add** ③ button.

The **Valco Valve Actuator Setup** dialog will appear.

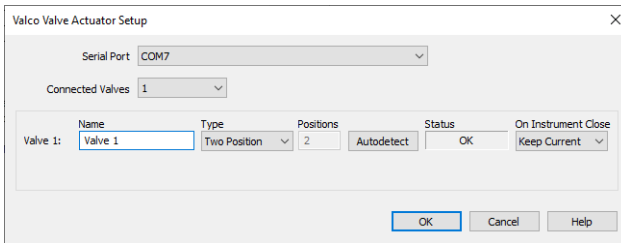

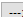


Fig 4: Valco Valve Actuator Setup

- Select the *Serial Port* and correct configuration of your valves.
- You might want to fill in custom valve names for each valve.

Note: The **Valco Valve Actuator Setup** dialog is more closely described in the chapter "**Valco Valve Actuator Setup**" on pg 10.

- The **ET, EH, EP, EHM, EMT, ECHM, ECMT** item ④ will appear in the *Setup Control Modules* list of the **System Configuration** dialog, with the number valves according the value set in the **Valco Valve Actuator Setup** dialog.
- Drag the individual Valve icons  from the *Setup Control Modules* ④ list on the left side of the **System Configuration** dialog to the desired *Instrument* ⑤ tab on the right side ⑥ (or use the  button ⑦ to do so).

Note: Individual Valves connected in the chain to the same Actuator can be configured on different Instruments.

- Set the *Ext. Start Dig. Input* and *Ready Dig. Output* numbers ⑧ for your acquisition card according to the wires being used for synchronization.

4 Using the Control Module

New **Valves** tabs (one for each valve assigned to the Instrument) are created in the **Method Setup** dialog. They can be accessed by switching to the desired valve in the **Select Valve** section on the top of the dialog.

4.1 Method Setup - Valves

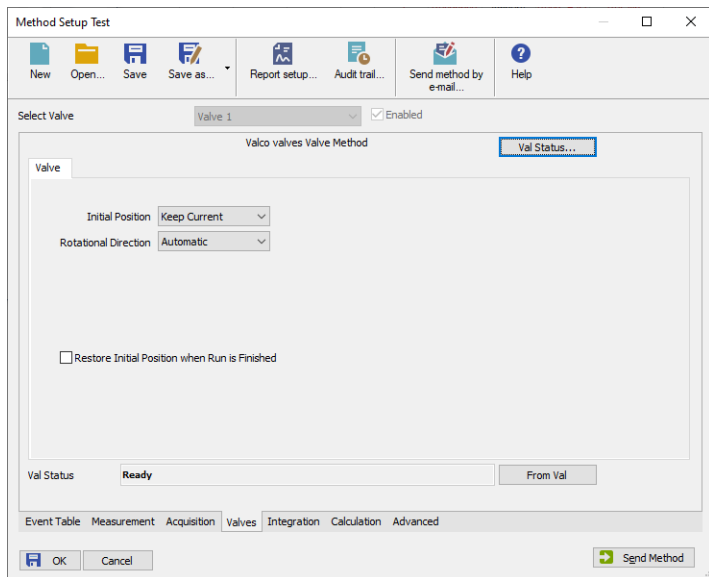


Fig 5: Method Setup - Valves

Initial Position

Specifies the initial position to which the valve is switched at the sending of the method. Possible values differ for the Two-position valves and Multiposition valves.

Possible values for the Two-position valves are: *Keep Current*, *A*, *B*.

Possible values for the Multiposition valves are: *Keep Current*, *1*, *2*, ..., *X*, where *X* is the number of positions set for the valve.

Keep Current option does not change the valve position at the sending of the method.

Rotational Direction

Sets the direction of valve position changes. The possible values are as follows: *Automatic*, *Forward*, *Reverse*. Setting the value to *Automatic* means that the valve switches from the current to the desired position the fastest way possible.

Note: This field is only present for Multiposition valves.

Restore Initial Position when Run is Finished

This checkbox specifies what will happen after the acquisition is finished. If checked, the valve will return to the state specified in the *Initial Position* field.

From Val

Acquires the status of the active valve from the device and sets the values obtained to the active valve's **Valves** tab of the **Method Setup** dialog.

Val Status

Invokes the **Hardware Configuration** dialog showing the COM port and list of valves connected to it via valve actuator. The firmware version of the actuator is shown next to each valve if the connection was successful, other possible values are *NOT CONNECTED* (the particular valve is not connected or switched on) and *ERROR* (error occurred during the valve test). The dialog is specific to the valve it is invoked from, e.g. the result shown will depend on the particular valve when invoked from the Instrument with two actuators configured on it.

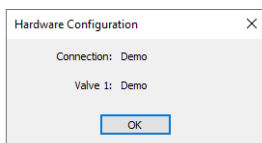


Fig 6: Hardware Configuration

4.2 Method Setup - Event Table

Event Table tab of the **Method Setup** dialog is the place from where the switching of valve positions is governed.

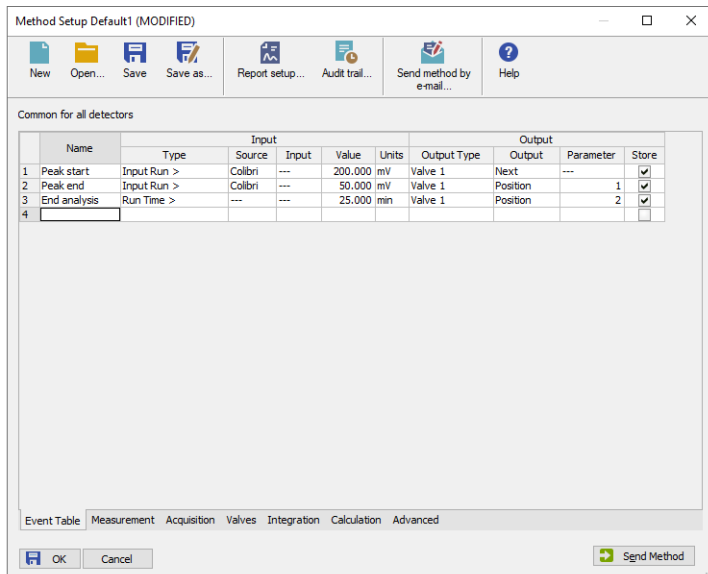



Fig 7: Method Setup - Event Table

The change of the valve position can be triggered by any kind of *Input* (analog signal absolute or relative value getting over or under the specified value, the analysis time passed, digital input, ...) and switched to any possible position. The valve to be switched is set in the *Output Type* field, while the *Output* field allows to select whether the valve should be switched to the next position (*Next*) or to any position defined by its number (*Position*). In that case, the *Parameter* field tooltip shows the possible range interval and the field itself allows to fill in the actual position number to switch into. Check the checkbox of the particular row in the *Store* column to log the event of switching the valve to the **Audit trail**.

4.3 Device Monitor

The window with the detector status can be invoked by the *Monitor - Device Monitor* command from the **Instrument** window or using the  *LC Monitor* icon. It displays the actual valve position for each valve and allows the user to change it if the analysis is not running.

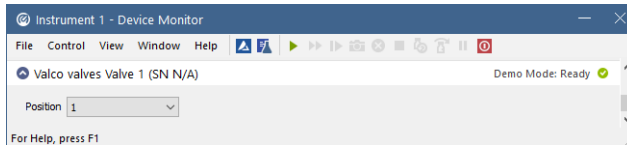


Fig 8: Device Monitor

Position

Shows the position of the valve and allows the user to set the desired valve position. During the run, this field is inactive (shaded).

4.4 Valco Valve Actuator Setup

The appearance of the **Valco Valve Actuator Setup** dialog depends on the number of connected valves set - the dialog has some functions different when more than one valve is selected.

Caution: If using more than one valve, do not connect them all at once. See the section concerning multiple connected valves.

One connected valve only

The screenshot shows the 'Valco Valve Actuator Setup' dialog box. At the top, there is a 'Serial Port' dropdown menu set to 'COM7' and a 'Connected Valves' dropdown menu set to '1'. Below these is a table with columns: Name, Type, Positions, Status, and On Instrument Close. The table contains one row for 'Valve 1' with the following values: Name: 'Valve 1', Type: 'Two Position', Positions: '2', Status: 'Autodetect', and On Instrument Close: 'Keep Current'. At the bottom of the dialog are three buttons: 'OK', 'Cancel', and 'Help'.

Name	Type	Positions	Status	On Instrument Close
Valve 1	Two Position	2	Autodetect	Keep Current

Fig 9: Valco Valve Actuator Setup - 1 connected valve

Serial Port

Sets the number of the serial port, to which the valves are connected.

Connected Valves

Sets the number of valves connected in chain to the actuator; up to 10 valves may be connected this way.

Name

Sets the name of the valve. This name will be used throughout the **Clarity** station.

Type

Select whether the controlled valve model is *Two Position* or *Multiposition*.

Positions

Sets the number of valve positions. The number can be only edited if the *Multiposition Type* was selected, otherwise the value is 2.

Autodetect

When invoked, checks the connection to the valve and displays the result in the *Status* field. Also tests the type of the valve (*Two Position* or *Multiposition*) and sets the correct type into the *Type* field.

Caution: User must ensure the Actuator Settings match the mounted valve to prevent its physical damage. The type and number of Valve positions are autodetected upon connection from the Valve Actuator. Those settings could be changed in **Clarity Valco Valve Actuator Setup**, but will not affect the connected Valve Actuator settings. Mismatch between configured and connected valve will be reported as communication error when opening **Clarity Instrument** window. It is necessary to fix this by connecting the valve with expected settings or change the **Clarity** configuration.

Status

Shows the current status of the particular valve. Possible states are:

OK - Valve is connected and ready.

NOT TESTED - Valve was not tested yet.

TESTING... - Valve is currently tested.

WRONG TYPE - Detected different valve type than it is selected.

NOT FOUND - Valve is not connected or is not switched on.

ERROR - Error occurred during valve test.

On Instrument Close

Defines the behavior of the valve when **Instrument** window is closed. The valve either does nothing (when *Keep Current* option is selected) or changes position to the selected one (other options).

Multiple connected valves

The screenshot shows the 'Valco Valve Actuator Setup' dialog box. At the top, 'Serial Port' is set to 'COM7' and 'Connected Valves' is set to '3'. Below this is a table with columns: Name, Type, Positions, Status, and On Instrument Close. There are three rows for Valve 1, Valve 2, and Valve 3. Each row has input fields for Name, Type, Positions, and Status, and a dropdown for On Instrument Close. A 'Test' button is located below the table. At the bottom of the dialog are 'OK', 'Cancel', and 'Help' buttons.

Name	Type	Positions	Status	On Instrument Close
Valve 1	Two Position	2	Identify	OK
Valve 2	Multiposition	16	Identify	OK
Valve 3	Multiposition	6	Identify	OK

If more valves are connected in chain, you have to identify each valve. Connect single valve and press "Identify" button of the particular valve.
Repeat identification for all valves one by one. Finally connect all valves in chain and press "Test" button.

Fig 10: Valco Valve Actuator Setup - more valves

The fields in this setup have the same meaning, except that the *Name*, *Type*, *Positions* and *Status* field are specific to the particular valve. The *Test* button is common to all connected valves and new *Identify* button is used.

Identify

Identifies the valve if more than valve is connected. A unique ID number is set to the valve by this procedure, which is consequently used in the

communication to address this particular valve. Also tests the type of the valve (*Two Position* or *Multiposition*) and sets the correct type into the *Type* field with the same functionality as the *Test* button.

Caution: When configuring more than one valve, it is necessary to connect and identify those valves one by one. Connect the first valve and press the *Identify* button in the first row. Then disconnect it, connect the second valve, press *Identify* button in the second row and so on. After all valves have been identified, it is possible to connect them all and test their status by the *Test* button.

Test

Functions in a similar way as the *Autodetect* button in single valve configuration - check connection to all identified valves and gets their current status.

5 Report Setup

The report of the valve part of the method can be enabled by checking the *Instrument Control* checkbox on the [Method](#) tab of the [Report Setup](#) dialog. Each valve configured on the Instrument has its own report section.

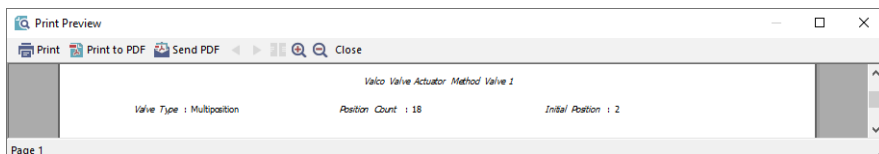


Fig 11: Report Setup

The type of the valve (multiposition vs. two-position), as well as the number of positions and the initial valve position are reported.

To report the actual events of the valve switches, it is necessary to check the checkboxes on the rows of the **Event Table** connected to the valves (so that these events are stored to the **Audit Trail**) and then check the *Chromatogram Audit Trail* option on the [Audit Trail](#) tab of the [Report Setup](#) dialog.

6 Troubleshooting

When the remedy for some problem cannot be discovered easily, the recording of communication between **Clarity** and the valves can significantly help the **DataApex** support to discover the cause of the problem.

The recording can be enabled by adding or amending the COMMDRV.INI file in the CFG subdirectory of the **Clarity** installation (C:\CLARITY\ by default). The file can be edited in any text editor (e.g. Notepad). Following section should be edited or added:

```
[COM1]
echo=on
textmode=on
filename=ViciValco_%D.txt
reset=off
```

Note: Instead of COM1 type the correct serial port used to communicate with the **VICI Valco Valves**. This port number is displayed when the *Val Status* button in the [Method Setup - Valve](#) or in the [Valco Valve Actuator Setup](#) dialog.

Note: %D (or %d) in the filename parameter means that the log will be created separately for each day. The *reset=off* parameter disables deleting the content of the log each time the station is started during the same day.

The created *.TXT files will greatly help in diagnosis of unrecognized errors and problems in communication.